


# DESIGNED

from the bottom up for your application

Ultra-low temperature freezer accessories

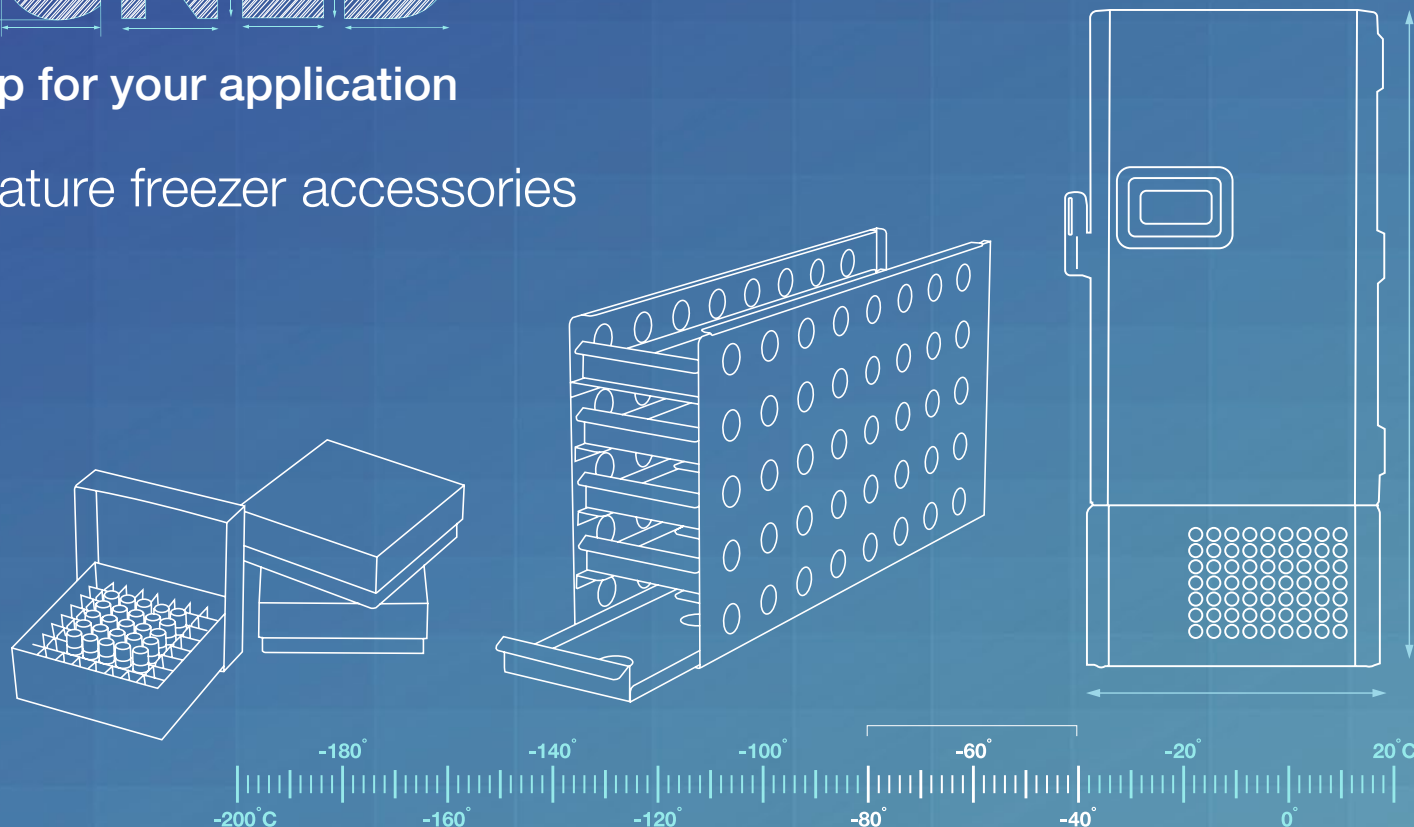

ULT accessories

## Racks and accessories guide for Thermo Scientific ULT freezers

Maximize storage space for ultimate  
sample protection

thermo  
scientific

Authorised Distributor



...řešení pro vaši laboratoř

TRIGON PLUS s.r.o.

T: +420 272 680 190, mail@trigonplus.cz, www.trigonplus.cz

Obchodní a servisní partner předních výrobců laboratorních zařízení

Zastoupení pro Českou a Slovenskou republiku

Akreditovaná kalibrační a zkušební laboratoř



thermo scientific

# Selecting the right accessories for your ULT freezer

Your Thermo Scientific™ ultra-low temperature (ULT) freezer is a serious investment in the long-term protection of your precious samples. It may not seem obvious at first, but the choice of your racks and accessories can help maximize your investment by optimizing operational efficiencies and reducing cost over the years.

## Which racks best suit your needs?

Selecting the appropriate racks for your application starts with two simple questions.

1. What kinds of samples are you storing?  The best choice of racks is determined by the nature of what you are storing.
2. Is your freezer an upright or chest unit?  The style of your freezer will determine the best selection of rack options. For instance, racks with side access and sliding drawers are available for upright freezers and offer unique advantages. The rack remains stationary and the drawers are light, making it easier to retrieve samples and minimize exposure to ambient conditions. Generally, racks are designed with specific dimensions and features, including handle placement, depending on the style of freezer. Understanding your storage needs up front and anticipating future requirements will help you determine the best choice of racks and accessories and help assure you'll gain maximum value from your ULT freezer.



# Rack and accessory options at a glance

Thermo Scientific™ ULT freezer racks and accessories are designed to integrate seamlessly with your existing equipment, enabling optimized performance and a more streamlined workflow. Discover tailored solutions that help to ensure your lab operates at peak efficiency.

- Precision-engineered for maximum storage and accessibility
- Corrosion-resistant materials to help ensure longevity and reliability
- Smooth surfaces protect both freezers and samples
- Versatile options for unique lab requirements

## Racks for boxes, microplates, and storage tubes



Adjustable side access rack



Chest freezer racks



Sliding drawer rack

Go to ordering table:

[Upright freezer racks](#) | [Chest freezer racks](#)

## Legacy and specialty racks



8-compartment container




Wire basket roller



10-compartment container

[Go to ordering table](#)

## Boxes and dividers

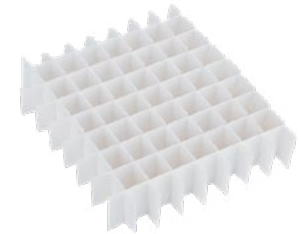

Fiberboard dividers



Fiberboard boxes


[Go to ordering table](#)

Various sizes and configurations available



## Contents

Racks for boxes and microplates	5
Racks for storage tubes	7
Legacy and specialty racks	9
Racking shelf kits	14
ULT chest freezer racks	15
Boxes and dividers	16

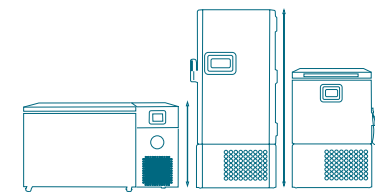

# Racks for boxes and microplates

## Suitable for 4-compartment ULT upright freezers


### Model series\* (4-compartment)

Description	Box type	Max box dimensions (L x W x H)	Height	Width	Depth	Storage capacities	Model series* (4-compartment)					Cat. No.
							TDE300	TSX400 TDE400	TSX500 TDE500	TSX600 TDE600	TSX700	
Front access sliding drawer rack	2 in./ 5.08 cm box	5.2 x 5.2 x 2.0 in.	11.9 in./30.2 cm	5.5 in./14 cm	25.79 in./65.5 cm	Boxes/Rack	25	25	25	25	25	920090
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	12	16	20	24	28	
						Boxes/Freezer	300	400	500	600	700	
Adjustable side access rack			11.6 in./29.5 cm	5.4 in./13.7 cm	26.75 in./67.9 cm	Boxes/Rack	25	25	25	25	25	1950520
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	12	16	20	24	28	
						Boxes/Freezer	300	400	500	600	700	
Front access sliding drawer rack	3 in./ 7.62 cm box	5.2 x 5.2 x 3.0 in.	11.9 in./30.2 cm	5.5 in./14 cm	25.79 in./65.5 cm	Boxes/Rack	15	15	15	15	15	920091
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	12	16	20	24	28	
						Boxes/Freezer	180	240	300	360	420	
Adjustable side access rack			11.6 in./29.5 cm	5.4 in./13.7 cm	26.75 in./67.9 cm	Boxes/Rack	15	15	15	15	15	1950521
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	12	16	20	24	28	
						Boxes/Freezer	180	240	300	360	420	
Front access sliding drawer rack	Microplates	5 x 5.3 x 0.6 in.	11.9 in./30.2 cm	5.5 in./14 cm	26.96 in./68.5 cm	Plates/Rack	35	35	35	35	35	1950642
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	12	16	20	24	28	
						Plates/Freezer	1,260	1,680	2,100	2,520	2,940	
Adjustable side access rack			11.9 in./30.2 cm	5.5 in./14 cm	25.5 in./64.8 cm	Plates/Rack	98	98	98	98	98	1950523
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	12	16	20	24	28	
						Plates/Freezer	1,176	1,568	1,960	2,352	2,744	
Adjustable side access rack	Microplates (standard or deep-well)	5 x 5.3 x 0.6 in. 5 x 5.3 x 1.8 in.	11.8 in./30.0 cm	5.5 in./14 cm	25.7 in./65.2 cm	Plates/Rack	147	147	147	147	147	1950592
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	12	16	20	24	28	
						Plates/Freezer	1,764	2,352	2,940	3,528	4,116	

Racks for boxes and microplates

[View imagery examples](#)

\* Also applies to the following series: 88000, 89000, FDE, TSU, and TSE.

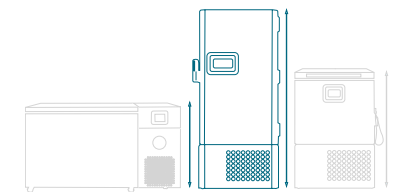

## Racks for boxes and microplates


### Suitable for 5-compartment ULT upright freezers

Description	Box type	Max box dimensions (L x W x H)	Height	Width	Depth	Storage capacities	Model series* (5-compartment)					Cat. No.
							TDE300	TSX400	TSX500	TSX600	TSX700	
								TDE400	TDE500	TDE600		
Front access sliding drawer rack	2 in./ 5.08 cm box	5.2 x 5.2 x 2.0 in.	9.5 in./24 cm	5.5 in./14 cm	26.9 in./68.3 cm	Boxes/Rack	20	20	20	20	20	398329
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	15	20	25	30	35	
						Boxes/Freezer	300	400	500	600	700	
Adjustable side access rack			9.4 in./23.9 cm	5.5 in./14 cm	26.75 in./67.9 cm	Boxes/Rack	20	20	20	20	20	1950518
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	15	20	25	30	35	
						Boxes/Freezer	300	400	500	600	700	
Front access sliding drawer rack	3 in./ 7.62 cm box	5.2 x 5.2 x 3.0 in.	9.5 in./24 cm	5.5 in./14 cm	26.9 in./68.3 cm	Boxes/Rack	15	15	15	15	15	398328
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	15	20	25	30	35	
						Boxes/Freezer	225	300	375	450	525	
Adjustable side access rack			9.4 in./23.9 cm	5.4 in./13.7 cm	26.75 in./67.9 cm	Boxes/Rack	15	15	15	15	15	1950519
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	15	20	25	30	35	
						Boxes/Freezer	225	300	375	450	525	
Front access sliding drawer rack	Microplates (standard or deep-well)	5 x 5.3 x 0.6 in. 5 x 5.3 x 1.8 in.	9.5 in./24 cm	5.5 in./14 cm	26.9 in./68.3 cm	Plates/Rack	28	28	28	28	28	1950639
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	15	20	25	30	35	
						Plates/Freezer	420	560	700	840	980	
Adjustable side access rack			9.4 in./23.9 cm	5.5 in./14.0 cm	25.5 in./64.8 cm	Boxes/Rack	77	77	77	77	77	1950522
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	15	20	25	30	35	
						Boxes/Freezer	1,155	1,540	1,925	2,310	2,695	

Racks for boxes and microplates

#### [View imagery examples](#)

\* Also applies to the following series: UxF, RLE, RDE, HFU-T, and HLE.

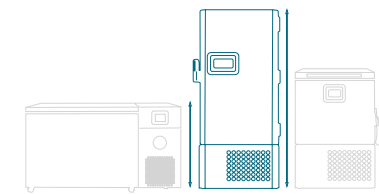


## Racks for storage tubes

Suitable for 4-compartment ULT upright freezers

Description	Tube type	Max box dimensions (L x W x H)	Height	Width	Depth	Storage capacities	Model series* (4-compartment)					Cat. No.
							TDE300	TSX400	TSX500	TSX600	TSX700	
								TDE400	TDE500	TDE600		
Front access sliding drawer rack	Matrix 200 µL, 2D or Matrix 500 µL, 2D	3.37 x 5.03 x 1.08 in.	11.9 in./30.2 cm	5.5 in./14.0 cm	26.96 in./68.5 cm	Tubes/Box	96	96	96	96	96	1950787
						Drawers/Rack	8	8	8	8	8	
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	12	16	20	24	28	
						Boxes/Freezer	672	896	1,120	1,344	1,568	
Front access sliding drawer rack	Matrix 750 µL, 2D	3.37 x 5.03 x 1.55 in.	11.9 in./30.2 cm	5.5 in./14.0 cm	26.96 in./68.5 cm	Tubes/Box	96	96	96	96	96	1950791
						Drawers/Rack	6	6	6	6	6	
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	12	16	20	24	28	
						Boxes/Freezer	504	672	840	1,008	1,176	
Front access sliding drawer rack	Matrix 1 mL, 2D	3.37 x 5.03 x 2.3 in.	11.9 in./30.2 cm	5.5 in./14.0 cm	26.96 in./68.5 cm	Tubes/Box	96	96	96	96	96	1950799
						Drawers/Rack	4	4	4	4	4	
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	12	16	20	24	28	
						Boxes/Freezer	336	448	560	672	784	
Front access sliding drawer rack	Matrix 1.4 mL, 1.8 mL, 2 mL, 2D	3.37 x 5.03 x 2.15 in.	11.9 in./30.2 cm	5.5 in./14.0 cm	26.96 in./68.5 cm	Tubes/Box	96	96	96	96	96	1950642
						Drawers/Rack	5	5	5	5	5	
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	12	16	20	24	28	
						Boxes/Freezer	420	560	700	840	980	
Front access sliding drawer rack	Nunc 5 mL, 2D	3.37 x 5.03 x 3.78 in.	11.9 in./30.2 cm	5.5 in./14.0 cm	26.96 in./68.5 cm	Tubes/Box	48	48	48	48	48	1950819
						Drawers/Rack	3	3	3	3	3	
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	12	16	20	24	28	
						Boxes/Freezer	252	336	420	504	588	

[View imagery examples](#)

\* Also applies to the following series: 88000, 89000, FDE, TSU, and TSE.

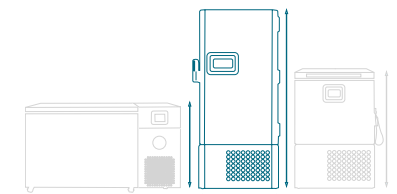

## Racks for storage tubes


Suitable for 5-compartment ULT upright freezers

Description	Tube type	Max box dimensions (L x W x H)	Height	Width	Depth	Storage capacities	Model series* (5-compartment)					Cat. No.
							TDE300	TSX400	TSX500	TSX600	TSX700	
								TDE400	TDE500	TDE600		
Front access sliding drawer rack	Matrix 200 µL, 2D or Matrix 500 µL, 2D	3.37 x 5.03 x 1.08 in.	9.5 in./24 cm	5.5 in./14.0 cm	26.96 in./68.5 cm	Tubes/Box	96	96	96	96	96	1950786
						Drawers/Rack	6	6	6	6	6	
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	15	20	25	30	35	
						Boxes/Freezer	630	840	1,050	1,260	1,470	
Front access sliding drawer rack	Matrix 750 µL, 2D	3.37 x 5.03 x 1.55 in.	9.5 in./24 cm	5.5 in./14.0 cm	26.96 in./68.5 cm	Tubes/Box	96	96	96	96	96	1950639
						Drawers/Rack	4	4	4	4	4	
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	15	20	25	30	35	
						Boxes/Freezer	420	560	700	840	980	
Front access sliding drawer rack	Matrix 1 mL, 2D	3.37 x 5.03 x 2.3 in.	9.5 in./24 cm	5.5 in./14.0 cm	26.96 in./68.5 cm	Tubes/Box	96	96	96	96	96	1950798
						Drawers/Rack	3	3	3	3	3	
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	15	20	25	30	35	
						Boxes/Freezer	315	420	525	630	735	
Front access sliding drawer rack	Matrix 1.4 mL, 1.8 mL, 2 mL, 2D	3.37 x 5.03 x 2.15 in.	9.5 in./24 cm	5.5 in./14.0 cm	26.96 in./68.5 cm	Tubes/Box	96	96	96	96	96	1950639
						Drawers/Rack	4	4	4	4	4	
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	15	20	25	30	35	
						Boxes/Freezer	420	560	700	840	980	
Front access sliding drawer rack	Matrix and Nunc 5 mL, 2D	3.37 x 5.03 x 3.78 in.	9.5 in./24 cm	5.5 in./14.0 cm	26.96 in./68.5 cm	Tubes/Box	24	24	24	24	24	1950818
						Drawers/Rack	2	2	2	2	2	
						Racks/Shelf	3	4	5	6	7	
						Racks/Freezer	15	20	25	30	35	
						Boxes/Freezer	420	560	700	840	980	

Racks for storage tubes

[View imagery examples](#)

\* Also applies to the following series: UxF, RLE, RDE, HFU-T, and HLE.

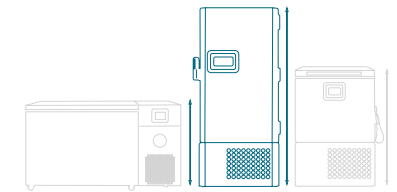

## Legacy racks

### Suitable for 4-compartment performance (legacy) ULT freezers


#### 4-compartment performance\* (legacy)

Description	Box type	Max box dimensions (L x W x H)	Height	Width	Depth	Storage capacities	902	905	906	907	Cat. No.	
							991	994	995			
Front access sliding drawer rack	2 in./5.08 cm box	5.2 x 5.2 x 2.0 in.	11.9 in./30.2 cm	5.5 in./14.0 cm	16.6 in./42.2 cm	Boxes/Rack	15				920094	
						Racks/Shelf	4					
						Racks/Freezer	16					
						Boxes/Freezer	240					
Adjustable side access rack			11.6 in./29.5 cm	5.3 in./13.5 cm	16.5 in./41.9 cm	Boxes/Rack	15				920015	
						Racks/Shelf	4					
						Racks/Freezer	16					
						Boxes/Freezer	240					
Front access sliding drawer rack			2 in./5.08 cm box	5.2 x 5.2 x 2.0 in.	11.9 in./30.2 cm	5.5 in./14.0 cm	21.97 in./ 55.8 cm	Boxes/Rack		20	20	920096
								Racks/Shelf		4	5	
								Racks/Freezer		16	20	
								Boxes/Freezer		320	400	
Adjustable side access rack	11.6 in./29.5 cm	5.3 in./13.5 cm			22.0 in./55.9 cm	Boxes/Rack		20	20	920018		
						Racks/Shelf		4	5			
						Racks/Freezer		16	20			
						Boxes/Freezer		320	400			
Adjustable side access rack	11.6 in./29.5 cm	5.3 in./13.5 cm			24.6 in./62.5 cm	Boxes/Rack				22	920022	
						Racks/Shelf				6		
						Racks/Freezer				24		
						Boxes/Freezer				528		
Front access sliding drawer rack	2 in./5.08 cm box	5.2 x 5.2 x 2.0 in.	11.9 in./30.2 cm	5.5 in./14.0 cm	26.3 in./66.8 cm	Boxes/Rack				25	1950726	
						Racks/Shelf				6		
						Racks/Freezer				24		
						Boxes/Freezer				600		

Legacy and specialty racks

[View imagery examples](#)

\* Also applies to Forma 900, Forma 700, TSE, and TSD.

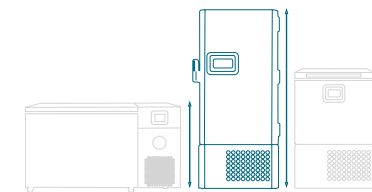

## Legacy racks

### Suitable for 4-compartment performance (legacy) ULT freezers


#### 4-compartment performance\* (legacy)

Description	Box type	Max box dimensions (L x W x H)	Height	Width	Depth	Storage capacities	902	905	906	907	Cat. No.				
							991	994	995						
Front access sliding drawer rack	3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.	11.9 in./30.2 cm	5.5 in./14.0 cm	16.6 in./42.2 cm	Boxes/Rack	9				920095				
						Racks/Shelf	4								
						Racks/Freezer	16								
						Boxes/Freezer	144								
Front access sliding drawer rack			3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.	11.9 in./30.2 cm	5.5 in./14.0 cm	22.0 in./55.9 cm	Boxes/Rack		12	12		920097		
								Racks/Shelf		4	5				
								Racks/Freezer		16	20				
								Boxes/Freezer		192	240				
Front access sliding drawer rack					3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.	11.9 in./30.2 cm	5.5 in./14.0 cm	26.3 in./66.8 cm	Boxes/Rack				15	1950727
										Racks/Shelf				6	
										Racks/Freezer				24	
										Boxes/Freezer				360	
Adjustable side access rack	3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.					11.6 in./29.5 cm	5.3 in./13.5 cm	16.5 in./41.9 cm	Boxes/Rack	9				920009
										Racks/Shelf	4				
										Racks/Freezer	16				
										Boxes/Freezer	144				
Adjustable side access rack			3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.			11.6 in./29.5 cm	5.3 in./13.5 cm	22.0 in./55.9 cm	Boxes/Rack		12	12		920017
										Racks/Shelf		4	5		
										Racks/Freezer		16	20		
										Boxes/Freezer		192	240		
Adjustable side access rack					3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.	11.6 in./29.5 cm	5.3 in./13.5 cm	24.6 in./62.5 cm	Boxes/Rack				15	1950521
										Racks/Shelf				6	
										Racks/Freezer				24	
										Boxes/Freezer				360	

Legacy and specialty racks

#### [View imagery examples](#)

\* Also applies to Forma 900, Forma 700, TSE, and TSD.

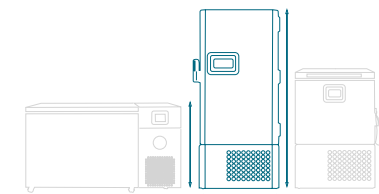

## Legacy racks

### Suitable for 5-compartment performance (legacy) ULT freezers


#### 5-compartment performance\* (legacy)

Description	Box type	Max box dimensions (L x W x H)	Height	Width	Depth	Storage capacities	ExF24086	ExF32086	ExF40086	ExF60086	Cat. No.				
Front access sliding drawer rack	2 in./5.08 cm box	5.2 x 5.2 x 2.0 in.	9.5 in./24.0 cm	5.5 in./14.0 cm	16.6 in./42.2 cm	Boxes/Rack	12				398332				
						Racks/Shelf	4								
						Racks/Freezer	20								
						Boxes/Freezer	240								
Adjustable side access rack			5.2 x 5.2 x 2.0 in.	5.2 x 5.2 x 2.0 in.	9.4 in./23.9 cm	5.5 in./14.0 cm	16.5 in./41.9 cm	Boxes/Rack	12				398324		
								Racks/Shelf	4						
								Racks/Freezer	20						
								Boxes/Freezer	240						
Front access sliding drawer rack					2 in./5.08 cm box	5.2 x 5.2 x 2.0 in.	9.5 in./24.0 cm	5.5 in./14.0 cm	22.0 in./55.9 cm	Boxes/Rack		16	16		398331
										Racks/Shelf		4	5		
										Racks/Freezer		20	25		
										Boxes/Freezer		320	400		
Adjustable side access rack	2 in./5.08 cm box	5.2 x 5.2 x 2.0 in.					9.4 in./23.9 cm	5.5 in./14.0 cm	22.1 in./56.1 cm	Boxes/Rack		16	16		398326
										Racks/Shelf		4	5		
										Racks/Freezer		20	25		
										Boxes/Freezer		320	400		
Front access sliding drawer rack			2 in./5.08 cm box	5.2 x 5.2 x 2.0 in.			9.5 in./24.0 cm	5.5 in./14.0 cm	26.3 in./66.8 cm	Boxes/Rack				20	1950728
										Racks/Shelf				6	
										Racks/Freezer				30	
										Boxes/Freezer				600	
Adjustable side access rack					2 in./5.08 cm box	5.2 x 5.2 x 2.0 in.	9.4 in./23.9 cm	5.5 in./14.0 cm	26.75 in./67.9 cm	Boxes/Rack				20	1950518
										Racks/Shelf				6	
										Racks/Freezer				30	
										Boxes/Freezer				600	

Legacy and specialty racks

#### [View imagery examples](#)

\* Also applies to Revco ExF and DxF and HERAFreeze HFU-B.

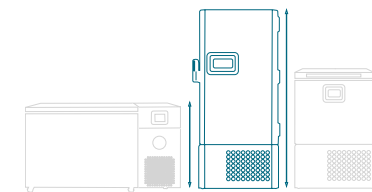

## Legacy racks


### Suitable for 5-compartment performance (legacy) ULT freezers

Description	Box type	Max box dimensions (L x W x H)	Height	Width	Depth	Storage capacities	5-compartment performance* (legacy)				Cat. No.				
							ExF24086	ExF32086	ExF40086	ExF60086					
Front access sliding drawer rack	3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.	9.5 in./24.0 cm	5.5 in./14.0 cm	17.4 in./44.2 cm	Boxes/Rack	9				920099				
						Racks/Shelf	4								
						Racks/Freezer	20								
						Boxes/Freezer	180								
Front access sliding drawer rack			3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.	9.5 in./24.0 cm	5.5 in./14.0 cm	22.0 in./55.9 cm	Boxes/Rack		12	12		398330		
								Racks/Shelf		4	5				
								Racks/Freezer		20	25				
								Boxes/Freezer		240	300				
Front access sliding drawer rack					3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.	9.5 in./24.0 cm	5.5 in./14.0 cm	26.3 in./66.8 cm	Boxes/Rack				15	1950729
										Racks/Shelf				6	
										Racks/Freezer				30	
										Boxes/Freezer				450	
Adjustable side access rack	3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.					9.4 in./23.9 cm	5.5 in./14.0 cm	16.1 in./40.9 cm	Boxes/Rack	9				398325
										Racks/Shelf	4				
										Racks/Freezer	20				
										Boxes/Freezer	180				
Adjustable side access rack			3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.			9.4 in./23.9 cm	5.5 in./14.0 cm	22.1 in./56.1 cm	Boxes/Rack		12	12		398327
										Racks/Shelf		4	5		
										Racks/Freezer		20	25		
										Boxes/Freezer		240	300		
Adjustable side access rack					3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.	9.4 in./23.9 cm	5.5 in./14.0 cm	26.75 in./67.9 cm	Boxes/Rack				15	1950519
										Racks/Shelf				6	
										Racks/Freezer				30	
										Boxes/Freezer				450	

Legacy and specialty racks

#### [View imagery examples](#)

\* Also applies to Revco ExF and DxF, and HERAFreeze HFU-B.

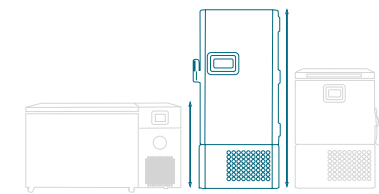


## Specialty racks

### For miscellaneous items

Description	Storage type	Height	Width	Depth	Cat. No.
8-compartment container	General freezer items	5.3 in./13.5 cm	7.0 in./17.8 cm	18.0 in./45.7 cm	830030
10-compartment container		5.3 in./13.5 cm	7.0 in./17.8 cm	22.6 in./57.4 cm	830032
Wire basket with roller for chest freezer*		11.8 in./29.9 cm	16.5 in./41.9 cm	12.3 in./31.2 cm	114048

\* Rail set required: Cat. No. 189001 for 13 cu. ft. chest freezers and Cat. No. 189000 for 20 cu. ft. chest freezers.

[View imagery examples](#)

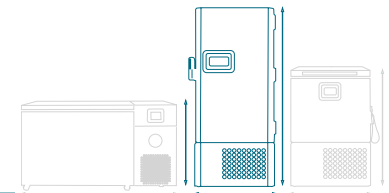


## Racking shelf kits

### For miscellaneous items

Description	Storage capacities	Model series and Cat. No.			
		TDE30086	TDE40086	TDE50086	TDE60086
		RSK300SD4	RSK400D4	RSK500SD4	RSK600SD4
Sliding drawer racks with boxes for one shelf (12.7 x 12.7 x 5.0 cm boxes with 100 count cell dividers)	Racks included	3	4	5	6
	Boxes included	75	100	125	150

Size-dependent options	Installation	TSX40086	TSX50086	TSX60086	TSX70086
Thermo Scientific™ Revco™ five inner door configuration	Factory	5D400	5D500	5D600	5D700
Shelf kit: includes one shelf and shelf support clips	Field	SK400TSX	SK500TSX	SK600TSX	SK700TSX

[View imagery examples](#)

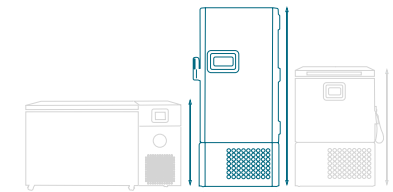


## ULT chest freezer racks

For miscellaneous items

Description	Box type	Max box dimensions (L x W x H)	Height	Width	Depth	Storage capacities	Model series			Cat. No.
							3 cu. ft.	13 cu. ft.	20 cu. ft.	
							TDEC06386	TDEC25286	TDEC39686	
Side access rack tower	2 in./5.08 cm box	5.2 x 5.2 x 2.0 in.	15.2 in./38.6 cm	5.5 in./14.0 cm	5.5 in./14.0 cm	Boxes/Rack	7			398184
						Racks/Freezer	9			
						Boxes/Freezer	63			
Side access rack tower	2 in./5.08 cm box	5.2 x 5.2 x 2.0 in.	26.5 in./67.3 cm	5.5 in./14.0 cm	5.5 in./14.0 cm	Boxes/Rack		12	12	398186
						Racks/Freezer		21	33	
						Boxes/Freezer		252	396	
Side access rack tower	3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.	16.1 in./40.9 cm	5.5 in./14.0 cm	5.6 in./14.2 cm	Boxes/Rack	5			398185
						Racks/Freezer	9			
						Boxes/Freezer	45			
Side access rack tower	3 in./7.62 cm box	5.2 x 5.2 x 3.0 in.	26.3 in./66.8 cm	5.5 in./14.0 cm	5.6 in./14.2 cm	Boxes/Rack		8	8	398187
						Racks/Freezer		21	33	
						Boxes/Freezer		168	264	
Side access rack tower	Microplates	5 x 5.3 x 0.6 in.	26.5 in./67.3 cm	5.5 in./14.0 cm	5.5 in./14.0 cm	Boxes/Rack		27	27	398188
						Racks/Freezer		21	33	
						Boxes/Freezer		567	891	

[View imagery examples](#)

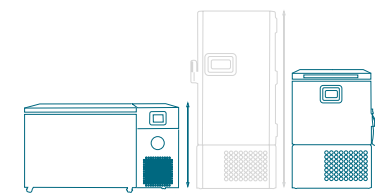

ULT chest freezer racks

## Boxes and dividers

### For ULT upright and chest freezers

Description	No. of dividers	Quantity	Color	Cat. No.
Thermo Scientific™ Nalgene™ polycarbonate 2 in. / 5.08 cm box for 1.2–2.0 mL vials	81	24/case	White	5026-0909
2 in. / 5.08 cm fiberboard box	0	12/pack	White	5954
2 in. / 5.08 cm fiberboard box	81	1	White	820002
2 in. / 5.08 cm fiberboard box	100	1	White	820109
3 in. / 7.62 cm fiberboard box	0	12/pack	White	5956
3 in. / 7.62 cm fiberboard box	81	1	White	820003
2 in. / 5.08 cm fiberboard box	0	12/pack	Multiple	1950878
2 in. / 5.08 cm fiberboard box	0	12/pack	Pink	1950879
2 in. / 5.08 cm fiberboard box	0	12/pack	Red	1950880
2 in. / 5.08 cm fiberboard box	0	12/pack	Yellow	1950881
2 in. / 5.08 cm fiberboard box	0	12/pack	Green	1950882
2 in. / 5.08 cm fiberboard box	0	12/pack	Blue	1950883
2 in. / 5.08 cm fiberboard box	0	12/pack	Black	1950884

#### Fiberboard dividers

Description	Quantity/pack	Cat. No.
Grid dividers, hold 100 x 12 mm tubes/vials	12	5958
Grid dividers, hold 81 x 13 mm tubes/vials	12	6212
Grid dividers, hold 64 x 14 mm tubes/vials	12	5960
Grid dividers, hold 49 x 16 mm tubes/vials	12	5959

[View imagery examples](#)

**thermo**  
scientific

Authorized Distributor



TRIGON PLUS s.r.o.

T: +420 272 680 190, mail@trigonplus.cz, www.trigonplus.cz

Obchodní a servisní partner předních výrobců laboratorních zařízení

Zastoupení pro Českou a Slovenskou republiku

Akreditovaná kalibrační a zkušební laboratoř

**thermo** scientific

Learn more at [thermofisher.com/ult](https://thermofisher.com/ult)

**For Laboratory Use. It is the customer's responsibility to ensure that the performance of the product is suitable for customers' specific uses or applications.** © 2024 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. **COL36239 0924**